

EQUIPMENT INSTALLATION

1

Position the base

2

Connect the Power Supply

3

Connect the heating device

4

Press UP button for 2 seconds to turn ON H401-T-PENNY

DEVICE SELECTION *

1

Press DOWN button for 2 seconds to enter in "Navigation Menu"

2

Press UP and DOWN buttons to enter in "Device Selection"

TYPE 000 DEVICE SELECTION

If the Type Number of your heating device is "Type 000", you have to manually input the control PID parameters (supplied by Okolab)

1

Press DOWN button for 2 seconds to enter in "Navigation Menu"

3

Read the heating device Type Number on the label

3

Press UP or DOWN button to select the heating device Type Number

2

Press UP and DOWN buttons to enter in "Device Selection"


3

Scroll the Type Number menu and press UP and DOWN buttons to select "Type 000"


*Remember to modify the Type Number if you change heating device.

TYPE 000 DEVICE SELECTION


4 Press UP and DOWN buttons to enter K1 subpages

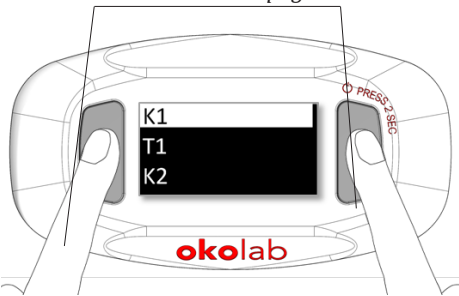

5 Press UP button to increase (or DOWN button to decrease) the K1 value

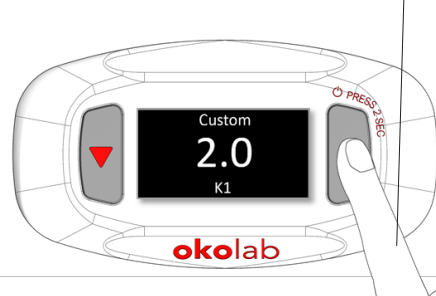


6 Press UP and DOWN buttons to confirm K1 value

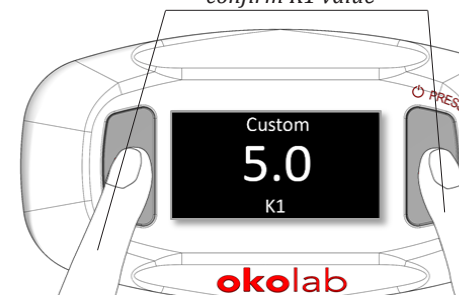


Repeat steps **4** **5** **6** to insert values for T1, K2, T2 and Delta T parameters

SET POINT TEMPERATURE


1 Press UP or DOWN button to select "Setpoint" from Homepage

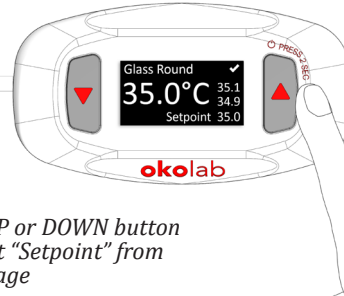

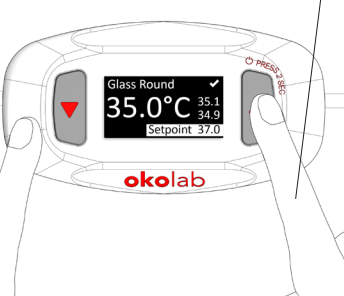

2 Press UP button to increase (or DOWN button to decrease) the Setpoint value

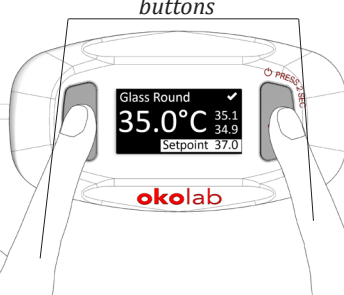

3 Confirm Setpoint value by pressing UP and DOWN buttons



As soon as the H401-T-PENNY stabilizes the temperature of the heating device, the Icon state changes from TRANSIENT  to NORMAL .